

Intelligent Type Peristaltic Pump

Model No. : LabV3-III

Flow Rate : 0.0053-1808mL/min



Single Channel



Dual Channel

Product Features

Industrial grade 4.3" true color LCD screen, touch screen control.

Mechanical keypad control pump start, stop, full speed and direction.

Dynamic display transferring status. Flow rate data, setting parameters and system configuration display in the same screen.

3 Kinds of working mode: fixed volume metering, fixed time and volume, timer start and stop, meet different transferring and dispensing request.

Intelligent calibration function and online micro adjusting function.

Screen Display

Real-time flow rate changeable by touching

Display area for setting parameter

Display area for dynamic working

Display area for running parameter

Function setting are changeable by touching

Fixed Volume Measurement Setting Interface

Dispensing Setting Interface

Time Start and Stop Setting Interface

Pump Head



EasyPump Series
(Pressure Adjustable)



EasyPump Series
(Fixed Pressure)



EasyPump Series
(Pressure Adjustable)



EasyPump Series
(Fixed Pressure)

Pump Head & Flow Rate

Pump Head	Tubing	ID×Wall thickness(mm)	mL / r	Flow Rate(mL/min) (LabV3-III, 0.1-350rpm)	Tubing Max.Pressure(Mpa)		Weight(kg)
					Intermittent	Continuous	
Single Channel EasyPump I/III	13#	0.8×1.6	0.053	0.0053~19	0.27	0.17	0.6
	14#	1.6×1.6	0.27	0.027~95			
	19#	2.4×1.6	0.55	0.055~193			
	16#	3.1×1.6	0.933	0.093~327	0.24	0.14	
	25#	4.8×1.6	1.967	0.197~688	0.14	0.10	
	17#	6.4×1.6	3.333	0.333~1167	0.10	0.07	
	18#	7.9×1.6	4.3	0.430~1505	0.27	0.17	
Single Channel EasyPump II/IV	15#	4.8×2.4	1.8	0.180~630			
	24#	6.4×2.4	2.733	0.273~957	0.24	0.14	
	35#	7.9×2.4	3.833	0.383~1342			
	36#	9.6×2.4	5.167	0.517~1808			
Dual Channel EasyPump V/VI	13#	0.8×1.6	0.0053	0.0053~19	0.27	0.17	
	14#	1.6×1.6	0.27	0.027~95			
	19#	2.4×1.6	0.55	0.055~193			
	16#	3.1×1.6	0.933	0.093~327	0.24	0.14	
	25#	4.8×1.6	1.967	0.197~688			

Pump Head	Tubing	Filling Volume(mL)	Filling Time(s)	Accuracy(±%)	Output(pcs/min)	Motor Speed(rpm)
EasyPump	13 [#]	0.4	1.2	0.8	27	377.36
	13 [#]	1	2.5	0.5	17	452.83
	14 [#]	2	1	0.5	30	444.44
	19 [#]	5	1.2	0.5	27	454.55
	16 [#]	7	1	0.5	30	450.16
	25 [#]	10	0.8	0.8	33	381.29
	25 [#]	15	1	0.5	30	457.55
	25 [#]	20	1.5	0.5	24	406.71
	17 [#]	30	1.2	0.5	27	450.05
	18 [#]	50	1.5	0.5	24	465.12
	15 [#]	15	1.2	0.5	27	416.67
	24 [#]	20	1.2	0.5	27	365.90
	35 [#]	30	1.2	0.5	27	391.34
	36 [#]	50	1.5	0.5	24	387.07
2*EasyPump	13 [#]	0.5	0.8	0.8	33	353.77
	13 [#]	1	1.2	0.5	27	471.70
	14 [#]	2	0.6	1	38	370.37
	14 [#]	5	1.5	0.5	24	370.37
	19 [#]	10	1.2	0.5	27	454.55
	16 [#]	15	1.2	0.5	27	401.93
	16 [#]	20	1.5	0.5	24	428.72
	25 [#]	30	1	0.5	30	457.55
	17 [#]	50	1	0.5	30	450.05
	18 [#]	70	1	0.5	30	488.37
	15 [#]	30	1.2	0.5	27	416.67
	24 [#]	50	1.2	0.5	27	457.37
	35 [#]	70	1.2	0.5	27	456.56
	36 [#]	100	1.5	0.5	24	387.07

Technical Specifications

Flow rate range	0.0053-1808mL/min
Flow rate resolution	0.01mL/min
Speed range	0.1~350rpm
Speed resolution	0.01rpm
Dispensing volume	0.01~9999mL
Volume resolution	0.01mL
Dispensing time	0.1~9999s
Pause time	0.1~9999s
Time resolution	0.01s
Repeat number	1-9999 times, setting "0" as infinite / continuous dispensing
Flow rate accuracy	<±0.5%
Back suction angle	0-360°
Motor type	Stepper motor
Display	4.3 inch-industrial grade-True color LCD screen
Control method	Touch screen and Mechanical keypad
Keypad lifetime	300,000 times
External control signal	0-5V,0-10V,4-20mA for option
Start/stop,change direction signa	Passive switch signal,such as:foot pedal switch,Active switch signal:5V-24V
Communication interface	RS232,RS485 support MODBUS protocol(RTU mode)
Output interface	Output motor working status(Open-Collector output)
Power supply	220V +10%50Hz/60Hz (standard),110V +10%50Hz/60Hz (optional)
Drive dimension (LxWxH)	259×157×237(L×W×H)
Drive Weight	4.40kg
Power consumption	<80W
condition temperature	0-40°C
Relative humidity	<80%

 Dimension Drawing (Unit:mm)

