

Model No. : OEM-Z216/EasyPump




Flow Data

Pump Head	Tubing	ID×Wall thickness(mm)	mL / r	Flow Rate(mL/min) (0.1-450rpm)	Tubing Max.Pressure(Mpa)		Weight(kg)
					Intermittent	Continuous	
Single Channel EasyPump I/III	13#	0.8×1.6	0.053	0.0053~23.85	0.27	0.17	0.6
	14#	1.6×1.6	0.27	0.027~121.5			
	19#	2.4×1.6	0.55	0.055~247.5			
	16#	3.1×1.6	0.933	0.093~419.85			
	25#	4.8×1.6	1.967	0.197~885.15	0.24	0.14	
	17#	6.4×1.6	3.333	0.333~1499.85	0.14	0.10	
	18#	7.9×1.6	4.3	0.430~1935	0.10	0.07	
Single Channel EasyPumpII/ IV	15#	4.8×2.4	1.8	0.180~810	0.27	0.17	
	24#	6.4×2.4	2.733	0.273~1229.85			
	35#	7.9×2.4	3.833	0.383~1724.85	0.24	0.14	
	36#	9.6×2.4	5.167	0.517~2325.15			
Dual Channel EasyPump V/VI	13#	0.8×1.6	0.0053	0.0053~23.85	0.27	0.17	
	14#	1.6×1.6	0.27	0.027~121.5			
	19#	2.4×1.6	0.55	0.055~247.5			
	16#	3.1×1.6	0.933	0.093~419.85			
	25#	4.8×1.6	1.967	0.197~885.15	0.24	0.14	

Motor Specification

Model No.	OEM-Z216/EasyPump
Motor type	DC geared motor
Power supply	DC24V
Speed	450rpm
Speed error	$\leq \pm 10\%$
Working environment	0-40°C, 80%RH

Dimension Drawing (Unit:mm)

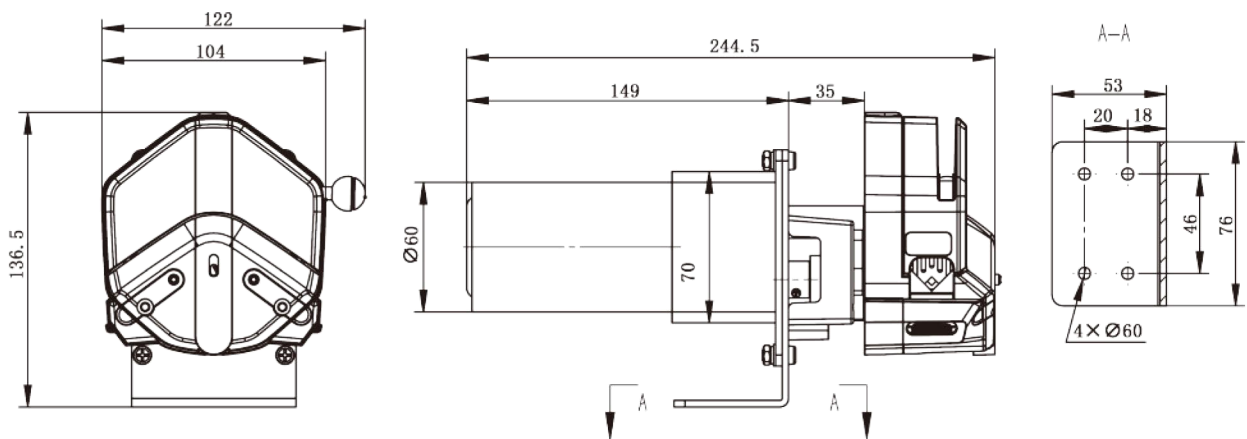

Tubing Installation Procedure




1. Turn the knob anticlockwise 180°, open the upper block.



2. Turn the rubbing wheel to adjust the lower tube clamp position, according to the tubing size.



3. Put the tube between rollers and upper block, tighten up the tubing lightly.

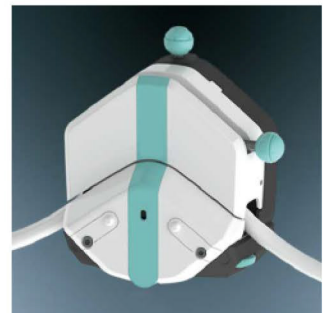

4. Turn the knob clockwise 180°, put the upper block back to original position, fasten the tube.